



Serrata
Pty. Ltd.

Tel: 02 9651 3033
Fax: 02 9651 2031
Email: serratalesley@ozemail.com.au
Web: www.serrata.com.au

Serrata Vacuum Pump

Instructions for Cat No 1014149



Read me

Vacuum pump 240V

Note: Oil containers are filled with light machine oil

Step by Step Instructions

1. Unpack the Vacuum Pump
2. Remove the oil bottles with care
3. Unscrew the white plug and remove
4. Fill the reservoir with oil using a funnel until the oil reaches the half way mark on the side window.
5. Replace and screw up the white plug
6. Remove the two black plugs
7. Untie the power lead
8. Plug the power plug into a 240V socket
9. Stay clear of the vacuum pump as a small amount of oil may be ejected the first time you turn it on
10. Turn on the switch
11. The two nozzles on the top of the Vacuum Pump are for:
 - a. Vacuum – Inlet (brass)
 - b. Exhaust – Outlet (black)
12. Turn off the Vacuum pump and store safely
13. Ensure the oil level is correct every time you use the Vacuum Pump

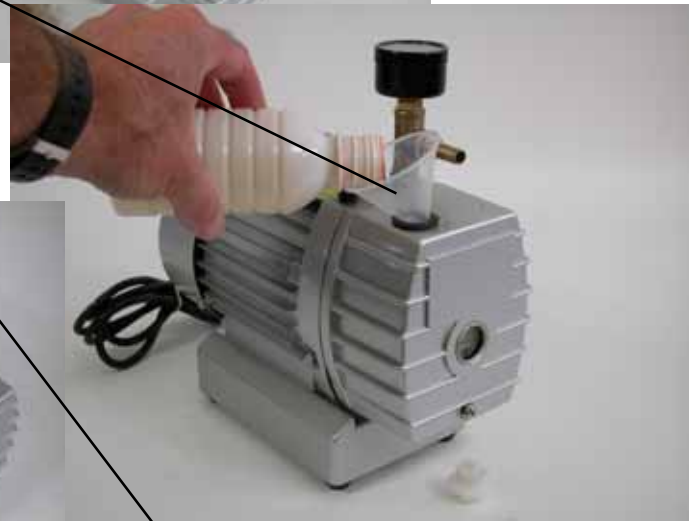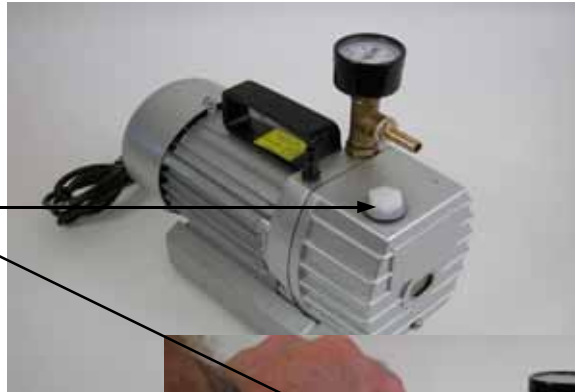

IMPORTANT. After evacuating a vessel and before switching off the pump, remove hose from inlet or otherwise release vacuum to prevent oil discharge.

Helping You Teach is Our Business!



IMPORTANT PLEASE NOTE:

This vacuum pump has been delivered to you pre-filled with oil.

This is because some schools were attempting to use the unit without reading the instructions, which stated that you needed to add oil before using it.

Before use, you only need to check that the oil level is above the red line but below the top of the sight glass.

The pump is also now fitted with a vacuum gauge on the inlet pipe, for greater accuracy and versatility.